


Precautions For Using Constant Climate Cabinet:

Mistake # 1: Preservation of medicine attached to internal condensing plate



The condensing plate is produced by the compressor. The cold source temperature is low. The medicine box will get wet slightly after contact, and it will freeze seriously. Keep a certain distance between the medicine box and the condensing plate to prevent contact.

Mistake # 2: Ceiling Fan Block

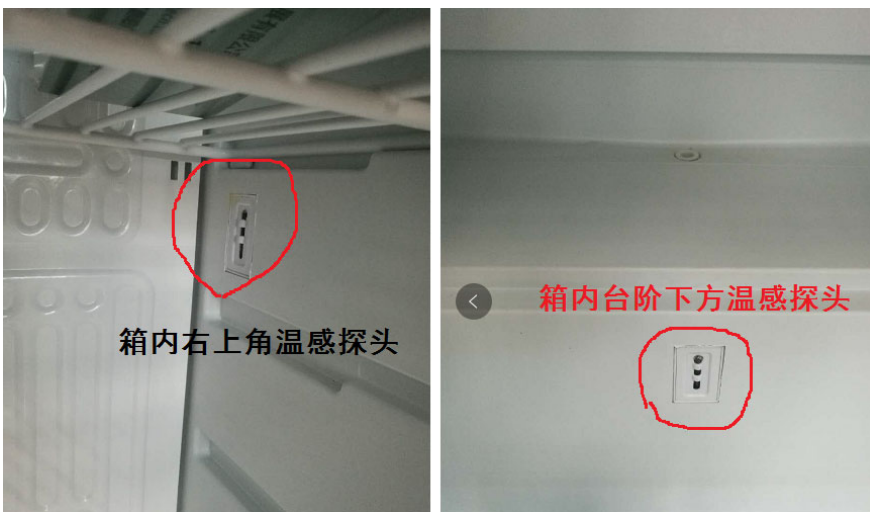

The ceiling fan generates circulating air to achieve the air circulation effect in the box, avoid over-temperature, and try to leave ventilation channels in the medicine placement.

Mistake # 3: Over-temperature or evaporator icing due to occlusion of temperature sensing probe.



Try to avoid the probe in the drug display to leave space for the probe, or due to the actual situation to extend the probe so as to achieve the correct temperature monitoring function.